

Combined SHIRPA and Dysmorphology TC

PLA_CSD_003

Purpose

SHIRPA and dysmorphology were originally always separate assessments. However they have recently been combined as assessments, so that they take place at the same time.

The purpose of the assessments is to examine mice for obvious physical characteristics, behaviors and morphological abnormalities.

Descriptions include abnormal locomotion/appearance/behavior/reflex reactions.

Experimental Design

- **Minimum number of animals** : 7M + 7F
- **Age at test**: Week 52
- **Sex**: We would expect the results of this test to show sexual dimorphism

Equipment

- Viewing Jar
- SHIRPA arena
- Grid above arena
- Click Box
- Geotaxis grid
- Tube for contact righting
- Photo and/or video camera

Procedure

1. Allow the mice to acclimatise to the phenotyping room for a period of 30 minutes prior to testing.
2. Throughout the test note any vocalisation, aggression, salivation or unexpected behaviours.
3. During this test, note any unexpected dysmorphological characteristics (morphological irregularities are recorded by video sequences or photos).
4. Place the mouse in a clear cylinder over a wire grid and observe for activity and tremors.
5. Transfer the mouse out of the cylinder by removing the metal plate/grid whilst positioning 30cm over an arena and record the transfer arousal.

6. Record the number of 10cm² squares the mouse moves into in the first 30 seconds in the arena (locomotor activity).
7. Allow the mouse to move freely around the arena whilst being observed for gait and tail elevation.
8. Hold the click box approximately 30cm above the arena and press the button, record the response of the mouse.
9. Pick up the mouse by the tail and observe for limp grasping and trunk curl. Trunk curl must only be recorded if the mouse curls forward without twisting its body, bending to one side is not scored as a trunk curl.
10. Place the mouse in a small transparent tube. Turn the tube quickly so the mouse is fully upside down and record if the mouse rights itself.
11. Record any vocalisation and/or aggression which were observed throughout the entire test

Notes

1. If wiping down with ethanol prior to the use of equipment, make sure no ethanol residue remains as the ethanol may affect the behaviour of the animals.
2. The validity of results obtained from behavioural phenotyping is largely dependent on methods of animal husbandry. It is important that individuals following this procedure are experienced and aware of the animal's welfare, and is familiar with the animal being tested, in order to reduce the anxiety levels of the animal prior to testing.
3. The majority of mouse behavioural studies are age/sex/strain dependent. It is important to keep these parameters comparable throughout a single experiment.
4. Environmental factors may contribute to the levels of anxiety within the mouse. The temperature, humidity, ventilation, noise intensity and light intensity must be maintained at levels appropriate for mice. It is essential that the mice be kept in a uniform environment before and after testing to avoid anomalous results being obtained.
5. It is recommended that all phenotyping experimentation is conducted at approximately the same time of day because physiological and biochemical parameters change throughout the day.
6. When a number of mice are tested continuously, residual odours from the equipment used in the preceding test may affect the test results. The floor and walls of the arena, ruler, and metal net should be wiped clean before introducing the next mouse. To prevent infection, the equipment should be washed with water at the completion of the day's tests. Some specific pathogen-free facilities use ultraviolet irradiation when tests are not being performed. Care needs to be taken, however, to ensure that ultraviolet irradiation does not crack any acrylate equipment covered with residual alcohol.

Parameters and Metadata

Hindlimb digit - size TCPLA_CSD_047_001 | v1.2

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both hindlimbs - increased size, One hindlimb - decreased size,
Both hindlimbs - fused, As expected, Both hindlimbs - decreased size,
One hindlimb - increased size, One hindlimb - fused,

Hindlimbs - size TCPLA_CSD_022_001 | v1.1

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: Both hindlimbs - decreased, One hindlimb only - increased,
Both hindlimbs - increased, One hindlimb only - decreased, As expected,

Head bobbing TCPLA_CSD_080_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: Absent, Present,

Transfer arousal TCPLA_CSD_031_001 | v1.2

simpleParameter

Req. Analysis: false **Req. Upload:** false **Is Annotated:** false

Options: As expected, Extended freeze, Immediate movement,

Teeth presence TCPLA_CSD_068_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: As expected, Not as expected,

Skin texture - ear TCPLA_CSD_064_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: Not as expected, As expected,

Skin texture - snout TCPLA_CSD_063_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: As expected, Not as expected,

Forelimb digit - number TCPLA_CSD_044_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both forelimbs - increased number, One forelimb - increased number, As expected, One forelimb - decreased number, Both forelimbs - decreased number,

Hindlimb digit - shape TCPLA_CSD_049_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both hindlimbs - not as expected, As expected, One hindlimb - not as expected,

Forelimb nail - length TCPLA_CSD_052_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: No data, One forelimb - decreased length, As expected, Both forelimbs - increased length, Both forelimbs - decreased length, One forelimb - increased length,

Skin texture - front paws TCPLA_CSD_065_001 | v1.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: Not as expected, As expected,

Size of squares in arena TCPLA_CSD_086_001 | v1.0

procedureMetadata

Req. Analysis: true Req. Upload: true Is Annotated: false

Unit Measured: cm^2

Skin color - snout TCPLA_CSD_057_001 | v1.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: As expected, Not as expected,

Limb grasp TCPLA_CSD_039_001 | v1.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: Absent, No data, Present,

Genitalia - morphology TCPLA_CSD_073_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected, No data,

Upper lip morphology TCPLA_CSD_075_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected,

Forelimb digit - shape TCPLA_CSD_048_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both forelimbs - not as expected, As expected, One forelimb - not as expected,

Coat - color - head TCPLA_CSD_006_001 | v1.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: No data, As expected, Not as expected,

Location of test TCPLA_CSD_082_001 | v1.1

procedureMetadata

Req. Analysis: false Req. Upload: true Is Annotated: false

Options: Open bench, LAF cabinet,

Skin color - tail TCPLA_CSD_061_001 | v1.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: Not as expected, No data, As expected,

Vocalization TCPLA_CSD_078_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Ears TCPLA_CSD_027_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Vibrissae - presence TCPLA_CSD_023_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Coat - hair texture / appearance - head TCPLA_CSD_015_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, No data, Not as expected,

Coat - hair distribution - head TCPLA_CSD_012_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: As expected, Not as expected,

Tail - length TCPLA_CSD_002_001 | v1.1

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: Long, As expected, No data, Short,

Hindpaw - size TCPLA_CSD_041_002 | v2.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: Both hindpaws - decreased, One hindpaw - decreased, As expected,
Both hindpaws - increased, One hindpaw - increased,

Coat - color - back TCPLA_CSD_005_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected, No data,

Hindlimbs - shape TCPLA_CSD_021_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: One hindlimb only - not as expected, Both hindlimbs - not as expected,
As expected,

Snout size TCPLA_CSD_028_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Decreased, Increased, As expected,

Lower teeth appearance TCPLA_CSD_070_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, No data, As expected,

Forelimb nail - number TCPLA_CSD_050_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both forelimbs - increased number, No data, Both forelimbs - decreased number, One forelimb - decreased number, As expected, One forelimb - increased number,

Skin color - back paws TCPLA_CSD_060_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Forelimbs - shape TCPLA_CSD_018_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both forelimbs - not as expected, As expected, One forelimb only - not as expected,

Contact righting TCPLA_CSD_077_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Present, Absent, No data,

Coat - hair texture / appearance - abdomen TCPLA_CSD_016_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected, No data,

Coat - color pattern - head TCPLA_CSD_009_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected, No data,

Unexpected behaviors TCPLA_CSD_037_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: Retropulsion, Jumping, None, Other, Circling,

Hindlimb nail - shape TCPLA_CSD_055_002 | v2.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: As expected, One hindlimb - not as expected, No data,
Both hindlimbs - not as expected,

Coat - color pattern - back TCPLA_CSD_008_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: No data, As expected, Not as expected,

Skin color - ear TCPLA_CSD_058_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected,

Head morphology TCPLA_CSD_026_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected,

Coat - color - abdomen TCPLA_CSD_007_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, No data, Not as expected,

Tail - presence TCPLA_CSD_001_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Absent, Present,

Forelimbs - size TCPLA_CSD_019_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both forelimbs - increased, One forelimb only - increased,
Both forelimbs - decreased, One forelimb only - decreased, As expected,

Trunk curl TCPLA_CSD_038_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Present, Absent, No data,

Tail - thickness TCPLA_CSD_003_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Decreased, Increased, As expected, No data,

Vibrissae - appearance TCPLA_CSD_024_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, No data, Not as expected,

Number of animals in cage TCPLA_CSD_083_001 | v1.2

procedureMetadata

Req. Analysis: false

Req. Upload: true

Is Annotated: false

Activity (body position) TCPLA_CSD_029_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Increased activity, Inactive, As expected,

Comments (in English) TCPLA_CSD_087_001 | v1.2

simpleParameter

Req. Analysis: false **Req. Upload:** false **Is Annotated:** false

Coat - hair distribution - abdomen TCPLA_CSD_013_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: As expected, Not as expected,

Aggression TCPLA_CSD_079_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: As expected, Excessive aggression,

Skin texture - tail TCPLA_CSD_067_001 | v1.0

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: As expected, Not as expected, No data,

Lower lip morphology TCPLA_CSD_076_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Coat - hair texture / appearance - back TCPLA_CSD_014_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: No data, Not as expected, As expected,

Hindlimb nail - length TCPLA_CSD_053_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both hindlimbs - decreased length, As expected, One hindlimb - decreased length, One hindlimb - increased length, Both hindlimbs - increased length, No data,

Forepaw - shape TCPLA_CSD_042_002 | v2.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: One forepaw - not as expected, As expected, Both forepaws - not as expected,

Startle response TCPLA_CSD_036_001 | v1.1

simpleParameter

Req. Analysis: false Req. Upload: false Is Annotated: true

Options: Present, No data, None,

Experimenter ID TCPLA_CSD_081_001 | v1.0

procedureMetadata

Req. Analysis: false Req. Upload: true Is Annotated: false

Images TCPLA_CSD_085_001 | v1.1

seriesMediaParameter

Req. Analysis: false Req. Upload: false Is Annotated: false

Tremor TCPLA_CSD_030_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Absent, Present,

Upper teeth appearance TCPLA_CSD_069_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, No data, As expected,

Hindlimb digit - number TCPLA_CSD_045_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: One hindlimb - increased number, Both hindlimbs - increased number,
Both hindlimbs - decreased number, As expected, One hindlimb - decreased number,

Forelimb nail - shape TCPLA_CSD_054_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: One forelimb - not as expected, As expected, Both forelimbs - not as expected,
No data,

Genitalia - presence TCPLA_CSD_071_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Absent, As expected,

Mouth morphology TCPLA_CSD_074_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Skin texture - back paws TCPLA_CSD_066_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected,

Hindlimb nail - number TCPLA_CSD_051_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both hindlimbs - decreased number, Both hindlimbs - increased number,
As expected, One hindlimb - decreased number, One hindlimb - increased number, No data,

Skin texture - whole body TCPLA_CSD_062_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Forelimbs - position TCPLA_CSD_017_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: One forelimb only - not as expected, As expected, Both forelimbs - not as expected,

Coat - color pattern - abdomen TCPLA_CSD_010_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: No data, As expected, Not as expected,

Hindlimbs - position TCPLA_CSD_020_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both hindlimbs - not as expected, As expected,
One hindlimb only - not as expected,

Locomotor activity TCPLA_CSD_032_001 | v1.2

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Unit Measured: Squares crossed

Head size TCPLA_CSD_025_001 | v1.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: Decreased, As expected, Increased,

Forepaw - size TCPLA_CSD_040_002 | v2.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: As expected, Both forepaws - decreased, One forepaw - decreased,
Both forepaws - increased, One forepaw - increased,

Gait TCPLA_CSD_033_001 | v1.0

simpleParameter

Req. Analysis: false Req. Upload: true Is Annotated: true

Options: Fluid movement, Lack of fluidity in movement,

Number of days since cage changed TCPLA_CSD_084_001 | v1.1

procedureMetadata

Req. Analysis: false

Req. Upload: false

Is Annotated: false

Tail - morphology TCPLA_CSD_004_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: No data, As expected, Not as expected,

Skin color - front paws TCPLA_CSD_059_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Forelimb digit - size TCPLA_CSD_046_001 | v1.1

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Both forelimbs - fused, As expected, One forelimb - fused,
One forelimb - increased size, Both forelimbs - decreased size,
One forelimb - decreased size, Both forelimbs - increased size,

Skin color - whole body TCPLA_CSD_056_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, Not as expected,

Coat - hair distribution - back TCPLA_CSD_011_001 | v1.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: Not as expected, As expected,

Hindpaw - shape TCPLA_CSD_043_002 | v2.0

simpleParameter

Req. Analysis: false

Req. Upload: true

Is Annotated: true

Options: As expected, One hindpaw - not as expected, Both hindpaws - not as expected,

Genitalia - size TCPLA_CSD_072_001 | v1.1

simpleParameter

Req. Analysis: false **Req. Upload:** true **Is Annotated:** true

Options: Not as expected, As expected, No data,

Tail elevation TCPLA_CSD_034_001 | v1.2

simpleParameter

Req. Analysis: false **Req. Upload:** false **Is Annotated:** false

Options: As expected, Dragging, No data, Straub / elevated tail,
